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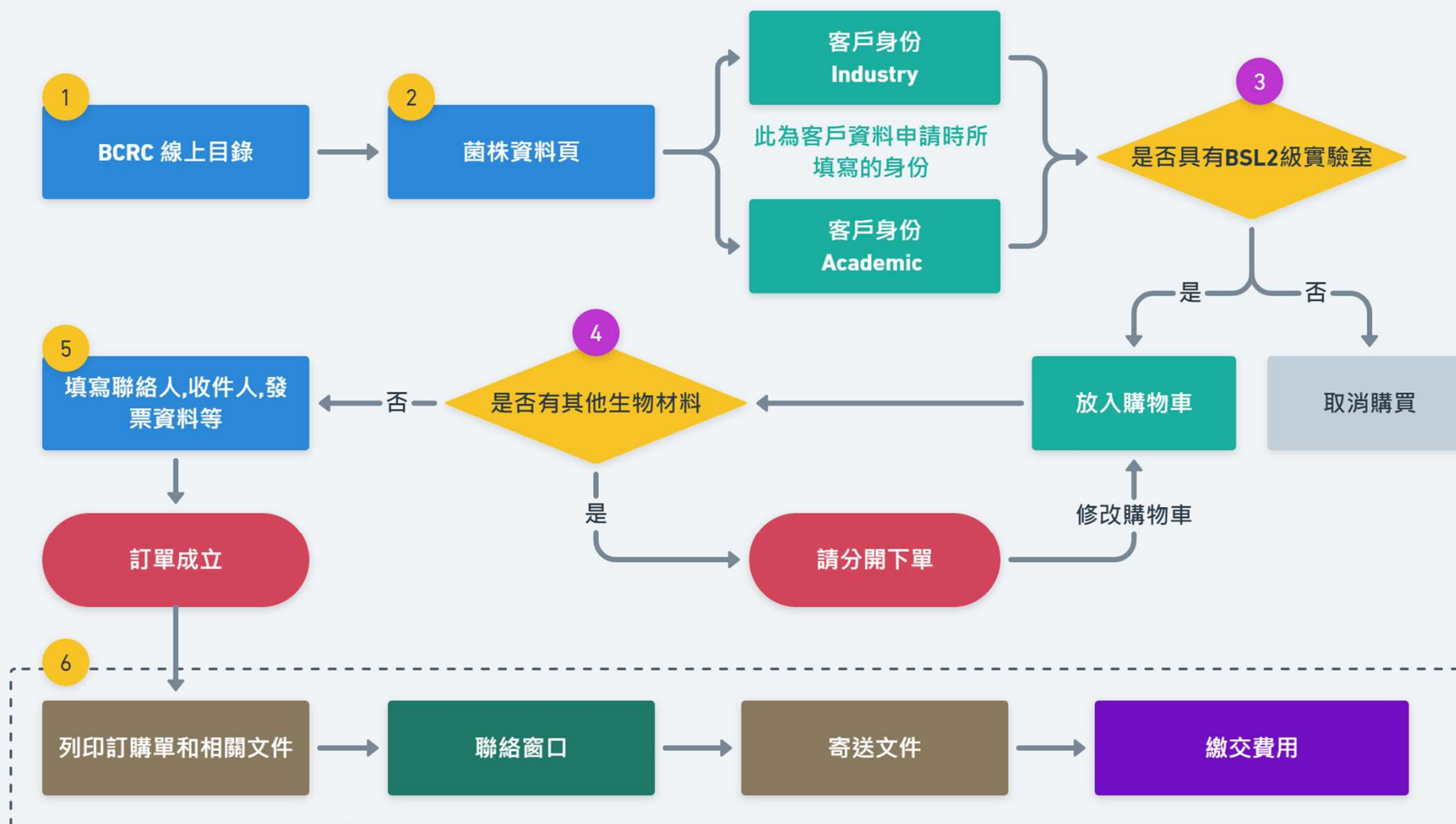
Home Reference Search Medium Search New Resources Reference Strain

Home

Search by Keyword

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Query

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Search by BCRC No.

ex.11633 21538 30506

More Search

Application strains

Lists for Applications

Anaerobe

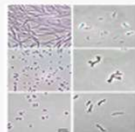
Nitrogen-fixing
Bacteria

Photosynthetic
Bacteria


Phosphorus-
solubilizing
Bacteria

Sulfur and Iron
Bacteria

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Bacteria
Search
A-Z

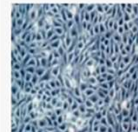Cell Line
Search
A-Z
iPS cell
List of discontinued cell
lines

Mushroom
Search
A-Z

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BCRC
Brain Collection Catalog
A Shipping Cart

Home Reference Search Medium Search New Resources Reference Strain

Home / Search in Cell Line / IPS cell / BCRC Content

[Return](#)

Cell Line

BCRC Number: SCB1034
Name: NTUH-IPSC-01-05
Type: Sendai virus reprogrammed human induced pluripotent stem cell
Species: Homo sapiens, Human
Age Stage: 45 years old
Gender: Female
Tissue: Peripheral blood mononuclear cell (PBMC)

Morphology: Colonies with smooth surfaces and well defined edges, ES-like
Split Ratio: Depend on culture vessel (10-20 colonies per 6cm dish)
Mycoplasma Test: Negative
Karyology: Normal 46, XX karyotypes

Incubation Temp(°C): 37°C
Incubation Condition: 5% CO₂ incubator
Culture Medium: 80% DMEM/F12 with 2mM L-glutamine adjusted to contain final sodium pyruvate + 1% NEAA + 1% GlutaMAX + 55uM 2-mercaptoethanol + 1% Penicillin-Streptomycin + 10ng/ml bFGF + 20% Knockout serum replacement (prepare the IPSC medium without bFGF, and then supplement with fresh bFGF when the medium is used)
Growth Property: Growth as colony-type
SubCulture Procedure: MEF based culture systems. FIRC recommends plating density for inactivated MEFs is 1x10⁵ cells per 10 cm dish (3.3x10⁵ cells 6 cm dish)

Days
Preparing MEF Culture Dishes
1. Cover the whole surface of each culture vessel with 0.1% gelatin and incubate the vessels for 30 minutes at 37°C or for 2 hours at room temperature.
2. Aspirate the gelatin solution from the gelatin coated culture vessel.
3. Wash the vessels with PBS.
4. Add complete IPSC Medium (pre-warm IPSC medium with b-FGF).
5. Use cell scraper cut the IPSCs to be picked into small pieces in a grid-like pattern.
6. Final transfer the IPSC pieces to the mouse embryonic fibroblasts (MEF) culture vessel containing IPSC medium.
7. Incubate the MEF culture plate containing the picked colonies in a 37°C incubator with a humidified atmosphere of 5% CO₂.
8. Change the medium every day.
9. DMEM/F12 + 25% Knockout serum replacement + 10% DMSO
Medium Renewal: Every day

Deposit History: National Taiwan University
Depositor: National Taiwan University
Deposit location: National Taiwan University
Category: IPS

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Characterization and validation to ensure the identity of the cell line: IPSC-01-05

Biosafety Level: 2
Use Restriction: After agreement of depositor, research use only
Sponsor: MOST-NCFB
Frozen Vial for academic NTD\$ 12000
Shipping/Field/Price: Frozen Vial for industry NTD\$ 36000

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